



# Thermo

ELECTRON CORPORATION  
Karl Fischer Titrator

Log # **KF1038**

## Overview

---

**Sample** PAPER-DISCS

**Industry** PAPER/PULP

**Group Number** 28

**Typical moisture** 1.7%

**Sample size** 1.0g

### Summary of procedure

### Method Parameters

<b>Instrument</b>	Turbo 2 Volumetric Karl Fische	<b>Titration Current</b>	
<b>Volumetric Reagent</b>	Reidel de Haen Composite 5	<b>Endpt Rate</b>	
<b>Volumetric Solvent</b>	MEOH	<b>Stirrer/blender speed</b>	3
<b>Pump Tubing</b>	SILICONE	<b>Extract time</b>	2.0min
<b>Vessel/Vessel Tubing</b>	STD 750mL	<b>Mode of analysis</b>	BKGD-SAMPLE-BKGD
<b>Endpt/Step Level</b>	8/5	<b>Vaporizer oven</b>	
<b>Endpt Time</b>	10sec	<b>Result units</b>	% wt/wt
<b>Anode Reagent</b>		<b>Source and date</b>	TW (031)
<b>Cathode Reagent</b>		<b>Cross-reference log</b>	KF1111

### Results

<b># of Trials</b>	5	<b>Mean Value</b>	1.66%	<b>C. V.</b>	4.8%
--------------------	---	-------------------	-------	--------------	------

### Comments