

## Thermo-Start™ PCR Kit

**Description:** Thermo-Start™ *Taq* DNA Polymerase is a chemically modified version of ThermoPrime *Taq* DNA Polymerase. It is completely inactive at room temperature, preventing the formation and subsequent amplification of non-specific products. The enzyme requires an **activation step at 95°C for 15 minutes**.

**Enzyme Source:** *Thermus aquaticus*

**Concentration:** 5 units/µl

**Unit Definition:** One unit of enzyme is defined as the amount that will incorporate 10nmoles of dNTPs into acid insoluble material in 30 minutes at 74°C under the analysis conditions below.

**Associated Activities:** Thermo-Start™ has 5' to 3' polymerization and exonuclease activity but lacks 3' to 5' exonuclease activity (proofreading).

**Storage Conditions:** Store Thermo-Start™ *Taq* DNA Polymerase at -20°C, in a constant temperature freezer for up to 12 months. Shipped on ice within the UK and on dry ice for international and within the US.

### Kit Contents:

Vial (cap color)	Pack Size	
	A	B
Thermo-Start <i>Taq</i> DNA Pol. (clear)	1 x 50µl	10 x 50µl
Thermo-Start PCR Buffer with MgCl <sub>2</sub> (purple)	1 x 1.25ml	10 x 1.25ml
dNTP Mix (brown)*	0.5ml	3 x 1.6ml

\*contains 5mM of each dNTP (dATP, dGTP, dCTP and dTTP)

**For Research Purposes Only**

1

**Example of Protocol:**

Mix and spin down the solutions prior to use

	Volume	Final Concentration 1X
Thermo-Start Taq DNA Pol. (5U/μl)	0.125μl	0.625 U
10X Thermo-Start PCR Buffer + 15mM MgCl <sub>2</sub>	2.5μl	1X (+1.5mM MgCl <sub>2</sub> )
dNTP Mix (20mM)	1μl	0.2mM of each nucleotide
Primer forward (10μM each)	1.25μl*	0.5μM*
Primer reverse (10μM each)	1.25μl*	0.5μM*
Water (PCR Grade)	variable	
DNA Template	0.5 - 10μl	0.5 - 125ng
Total Volume	25 μl	

\*Scale up or down the volume and concentration as appropriate

Note: These recommendations are intended as basic guidelines.

**Example of Program:**

	Temp.	Time	Number of cycle
<b>Initial Denaturation</b>	<b>95°C</b>	<b>15 min</b>	<b>1 cycle</b>
Denaturation	95°C	20 sec	30 to 40 cycles
Annealing	50-65°C	30 sec	
Extension**	72°C	60 sec	
Final Extension	72°C	5 min	1 cycle

\*\*Increase length of time in proportion to size of amplicon, Taq DNA Polymerase extends at approximately 1000 bp/min.

**Incremental Activation:**

For extra stringency, the enzyme can be activated gradually during the PCR in a series of steps. The initial activation step is replaced by longer (2 minutes) denaturation steps for the first 7–8 cycles of the reaction.

<b>Analysis Conditions:</b>	25mM	TAPS, pH 9.3 (at 25°C)
		[tris-(hydroxymethyl)-methyl-amino-propane sulfonic acid, sodium salt]
	50mM	KCl
	2mM	MgCl <sub>2</sub>
	1mM	β-mercaptoethanol
	200µM	of each: dATP, dGTP, dTTP
	100µM	[α- <sup>32</sup> P]-dCTP (0.05 to 0.1 Ci/mmol)
	1.25µg/µl	activated salmon sperm DNA

Water added to a total volume of 50µl. Incubated at 74°C for 10 minutes. The enzyme is first treated with a 15 minute activation step at 95°C. The amount of incorporated dNTPs is determined by trichloroacetic acid precipitation.

<b>Storage Buffer:</b>	100mM	KCl
	20mM	Tris-HCl, pH 9.2 (at 25°C)
	0.1mM	EDTA (ethylenediaminetetraacetic acid)
	1mM	DTT (dithiothreitol)
	0.5%	Tween <sup>®</sup> 20
	0.5%	Nonidet <sup>®</sup> P40
	50% (v/v)	Glycerol

<b>Ordering Information:</b>	AB-1908/A/N	Thermo-Start PCR Kit	250 units
	AB-1908/B/N	Thermo-Start PCR Kit	2500 units

All sizes are supplied with 10X Reaction Buffer IV containing 15mM MgCl<sub>2</sub> and a 20mM dNTP Mix.

## Troubleshooting

For troubleshooting, see [www.abgene.com/troubleshoot.asp](http://www.abgene.com/troubleshoot.asp) or contact Thermo Fisher Scientific (ABgene) TechSupport at [abgene.techsupport@thermofisher.com](mailto:abgene.techsupport@thermofisher.com)

UK TechSupport, call +44 (0) 1372 840 410

**For all other regions, please contact your local Thermo Fisher Scientific (ABgene) office / distributor.**

Use of this product is covered by one or more of the following US patents and corresponding patent claims outside the US: 5,079,352, 5,789,224, 5,618,711, 6,127,155, 5,677,152, 5,773,258, and claims outside the US corresponding to US Patent No. 4,889,818. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. No right under any other patent claim (such as the patented 5' Nuclease Process claims in US Patents Nos. 5,210,015 and 5,487,972) and no right to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is conveyed expressly, by implication, or by estoppel. This product is for research use only. Diagnostic uses under Roche patents require a separate license from Roche. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

**For Research Purposes Only**

3