

# Installation Request Checklist

## Surveyor MSQ

1. To insure proper site requirements for installation, please complete this checklist.  
Refer to the **Pre-Installation guide** for more information about each step.

- All laboratory remodeling has been completed.
  - Your Surveyor MSQ system is on site.
  - The entrance to the laboratory is at least 70 cm (28") wide.
  - The principal operator will be available during the installation period.
  - The workbench can support the **load** of the system (100 kg / 268 lbs) and is free from vibration.
  - The laboratory can be maintained at a constant temperature, between 15° and 35°C (63° and 95°F).
  - The relative humidity is between 40% and 80%, with no condensation.
  - The system work area is free from magnetic disruption and electrostatic discharge.
  - There is a telephone near the system.
  - There is a suitable **exhaust** system.
  - All relevant safety regulations are in compliance.
  - HPLC is new or recently cleaned, producing pulse free flow from 100-2000 micro-liters per minute.
  - Solvents:** HPLC grade water, methanol, acetonitrile and ammonium hydroxide are available.
  - Sufficient bench **space** is available for all the equipment, please list width: \_\_\_\_\_ depth: \_\_\_\_\_
  - The **power** outlets have been measured. Please note voltages: \_\_\_\_\_ volts.
  - The **power** outlets are the correct configuration. Please note NEMA types: \_\_\_\_\_
  - Nitrogen **gas** supply regulated at 75 psig (capable of 700 l/hr flow) with a purity of <sup>3</sup> 99% is available.
- Have any special acceptance specifications been agreed within the contract? Yes  No   
If **YES**, please attach full details of specifications.
- Is there any additional equipment that needs to be interfaced to the system? Yes  No   
If **YES**, please supply details.

## 2. Request Installation

To **request installation**, please review this form for accuracy and **mail or fax** to your local sales/service office:  
Schaumburg, Illinois: **847-310-0498**, San Jose, California: **408-965-6123**, Piscataway, New Jersey: **732-981-0029**,  
Woodstock, Georgia: **770-516-6916**

Name \_\_\_\_\_

Company \_\_\_\_\_

Address \_\_\_\_\_

City/State/Zip \_\_\_\_\_

Phone \_\_\_\_\_

Please note: We reserve the right to invoice against the engineers' time if the installation requirements are not met on date of install.

Signature: \_\_\_\_\_

Date: \_\_\_\_\_

**Thermo**Finnigan