

## Installation Request Checklist Trace DSQ

**1. To insure proper site requirements for installation, please complete this checklist.  
Refer to the Pre-Installation guide for more information about each step.**

- All laboratory remodeling has been completed and your Trace DSQ is on site.
- The entrance to the laboratory is at least 102cm (40") wide.
- The principal operator will be available during the installation period.
- The workbench is at least 2 m (6 ft) across and can support the load of the system (262 lbs or 119 kg) and is free from vibration.
- The laboratory can be maintained at a constant temperature, between 15° and 31°C (59° and 88°F).
- The relative humidity is between 40% and 80%, with no condensation.
- The system work area is free from magnetic disruption, electrostatic discharge and is properly lit.
- There is a telephone near the system.
- There is a suitable exhaust system.
- All relevant safety regulations are in compliance.
- The recommended purity solvents are available.
- Sufficient bench space is available for all the equipment, please list width: \_\_\_\_\_ depth: \_\_\_\_\_
- The power outlets have been measured. Please note voltages: \_\_\_\_\_ volts.
- The power outlets are the correct configuration. Please note NEMA types: \_\_\_\_\_
- The following gases are on-site with gas lines installed and appropriate regulators available.  
Please list gases and purity: \_\_\_\_\_

Have any special acceptance specifications been agreed within the contract? Yes  No   
If YES, please attach full details of specifications.

Is there any additional equipment that needs to be interfaced to the system? Yes  No   
If YES, please supply details.

### 2. Request Installation

To request installation, please review this form for accuracy and mail or fax to your local sales/service office:  
Schaumburg, Illinois: 847-310-0498, San Jose, California: 408-965-6123, Piscataway, New Jersey: 732-627-0260,  
Woodstock, Georgia: 770-516-6916

Name \_\_\_\_\_

Company \_\_\_\_\_

Address \_\_\_\_\_

City/State/Zip \_\_\_\_\_

Phone/Fax \_\_\_\_\_

Please note: We reserve the right to invoice against the engineers' time if the installation requirements are not met on date of install.

Signature: \_\_\_\_\_

Date: \_\_\_\_\_

Email \_\_\_\_\_