Instructions for Safe Use of IVD CE marked Finnpiquette® and Finntip®

The CE mark covers the combination (1514141, Finnpiquette -Finntip Compatibility Table) of CE marked Finnpiquette and the CE marked Finntip. Any other pipette - tip combination has to be validated by the user. In IVD use, the Finnpiquette is to be used by **trained personnel**, who have adequate laboratory skills. The pipette shall be **regularly serviced** according to the instructions for use.

Intended Use: The intended use of the Finnpiquette (type IVD) is to transfer liquids.

Performance: When leaving the factory, the pipette is tested against the **factory specifications**. The performance in normal use may vary (due to; temperature, pressure, humidity, operator, liquid, type of tip etc.). The user may also check the performance of the pipette with water against **limits of the standard ISO 8655** (See Instructions for Use). The user shall determine the **performance and applicability** of the pipette and tip combination **for the IVD application** and determine the required interval for performance checking. If the pipetting performance is critical to the medical diagnosis the diagnostic test result has to be ensured with an alternative test.

Prior to use Read the Instructions for Use. The performance of each new pipette needs to be checked and documented prior to use. The performance of each pipette after service needs to be checked and documented prior to use. Check that the package of sterile tips is unbroken (to ensure sterility). Check that the tip is complete and unbroken. Check that the filtertips have filters in place and the filters are unbroken or not tilted. Do not reuse the tips. Before use, check that there is a tip attached to the pipette.

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